

NIPRO CELL GATCH

For more information : http://www.funakoshi.co.jp/exports_contents/80659

NIPRO CELL GATCH is a protective pouch for cryotubes, made of special film, which is durable even in liquid nitrogen. During storage of cells in liquid nitrogen or when thawing cells, damage to the cryotubes is one of the biggest issues as this causes contamination of liquid nitrogen. NIPRO CELL GATCH prevents such contamination or sample scattering.

Features

- Special film, durable in liquid nitrogen
- Transparent; can see inside the cryotube
- Easy handling




Handling Procedure

■ Freezing

- (1) Put a cryotube into a pouch.



- (2) Seal two points within two lines of a pouch by using impulse sealer. Be careful not to wrinkle the sealing side.

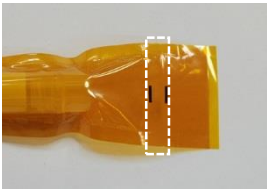

- (3) Put the pouches into a cane carefully not to damage. Using a cryo-sleeve is recommended to avoid drop of tubes from cane.



■ Thawing

- (1) Take a cane from liquid nitrogen and remove pouch carefully.
- (2) Soak the whole pouch in a water bath for thawing. After thawing, cut the pouch with scissors and take cryotube out.

Product Information

[Manufacturer : KSK]

Product Name	Size	Code	Storage
NIPRO CELL GATCH	20 sheets	87-449	RT

2016

Disposable Hemocytometer

For more information : http://www.funakoshi.co.jp/exports_contents/80019


made in Japan



High quality and Cost effective



- Disposable cell counting chamber slide
- Designed for rapid loading
- No washing, no labor, save time
- Same grid as standard hemocytometer
- High calculation accuracy
- Made of durable plastic

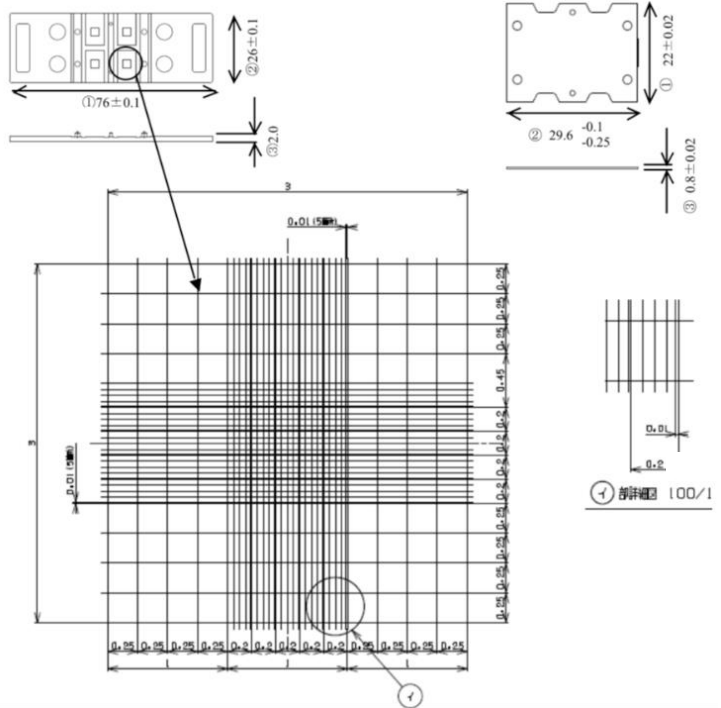

Free samples available !

Product name	Code	Unit
Disposable Hemocytometer (4-chambers)	521-10	50 pieces / 1 case
		200 pieces / 4 cases

NOTE

※ All products here are research use only, not for diagnostic use.
※ Specs might be changed for improvement without notice.

※ Company name and product name are trademark or registered mark.
※ Please contact your local distributors for orders, quote request and inquiry.

Your Local Distributor

Biolinkk

Building NO: 1333, 4th Floor, Office No:412,
Durga Chamber, Desh Bandhu Gupta Road,
Karol Bagh, New Delhi- 110005, India
Phone: +91-11-45035753
email: customerservice@biolinkk.com



funakoshi Co., Ltd.

Address: 9-7 Hongo 2-Chome, Bunkyo-ku,
Tokyo 113-0033 JAPAN
Phone : +81-3-5684-6296
Fax : +81-3-5684-6297
Email : export@funakoshi.co.jp