
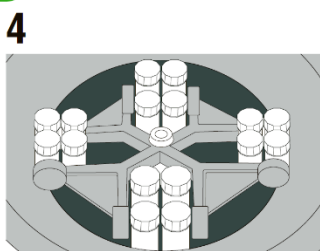
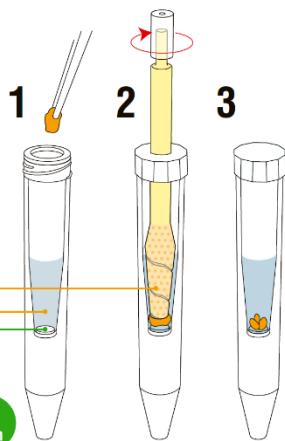


Convenient and Safe! Disposable Homogenizer BioMasher V

For more information : http://www.funakoshi.co.jp/exports_contents/80608

BioMasher V is a convenient disposable homogenizer. This is suitable for extraction of DNA, RNA and protein from samples.

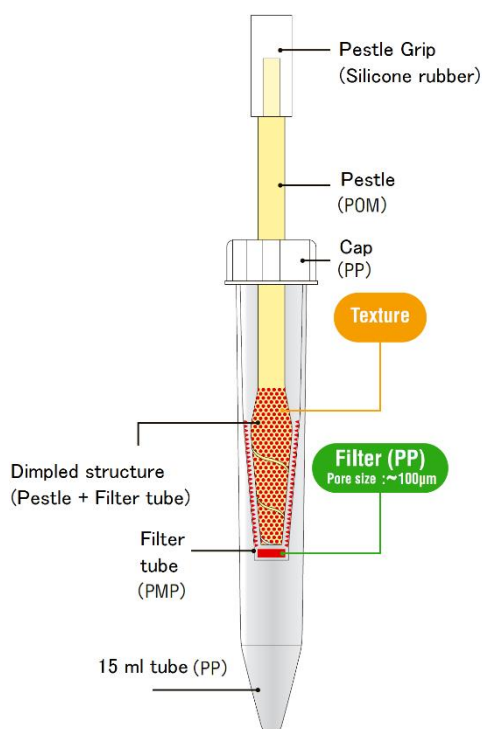

Features

- High homogenization efficiency
 - By texture on both pestle and inside of filter tube.
- High collection amount
 - Enlarged type of BioMasher III.
- Heat can be controlled
- Centrifuge required
- PP : Autoclavable / POM : non-autoclavable
- EOG-Sterilized or non-sterilized product are available.
- Much higher homogenization efficiency by using PowerMasher II (Pestle Motor).

Handling procedure

1. Put samples into tube and insert pestle.
2. Grinding samples by pestle by hand or PowerMasher II.
3. Remove pestle.
4. Centrifuge at 6,000 to 10,000 x g for 10 - 30 seconds.

PowerMasher II

PowerMasher II is a pestle motor, optimized for BioMasher II, III, V and SP series.

Features

- Increase homogenization efficiency dramatically.
- Works on two AA batteries (Accessories).



Check a movie at
http://www.funakoshi.co.jp/exports_contents/80608

Product Information

■ BioMasher V

[Manufacturer : NIP]

Product Name	Size	Code
BioMasher V (non-sterilized)	20 sets	891390
	100 sets	891391
BioMasher V (EOG-sterilized)	20 sets	891392
	100 sets	891393

■ PowerMasher II (Pestle motor)

[Manufacturer : NIP]

Product Name	Size	Code
PowerMasher II	1 piece	891300

NOTE

- * All products on here are for research use only, not for diagnostic use.
- * Products might be changed for improvement without notice.

- * All brand names and product names are registered trademarks of their respective companies.

- * Please contact your local distributors for orders, quote request and any inquiry.

Your Local Distributor

Biolinkk

Building NO: 1333, 4th Floor, Office No:412,
 Durga Chamber, Desh Bandhu Gupta Road,
 Karol Bagh, New Delhi- 110005, India
 Phone: +91-11-45035753
 email: customerservice@biolinkk.com



funakoshi Co., Ltd.

Address: 9-7 Hongo 2-Chome, Bunkyo-ku,
 Tokyo 113-0033, JAPAN
 Phone: +81-3-5684-6296
 Fax: +81-3-5684-6297
 Email: export@funakoshi.co.jp