

Easy to Use! Just Stick it on your Slide



PAP Barrier Seal for IHC

Only One


For more information : https://www.funakoshi.co.jp/exports_contents/520060?FL



Having troubles with PAP pen enclosure?
Now you are free from peeling off, drying time or liquid leakage!

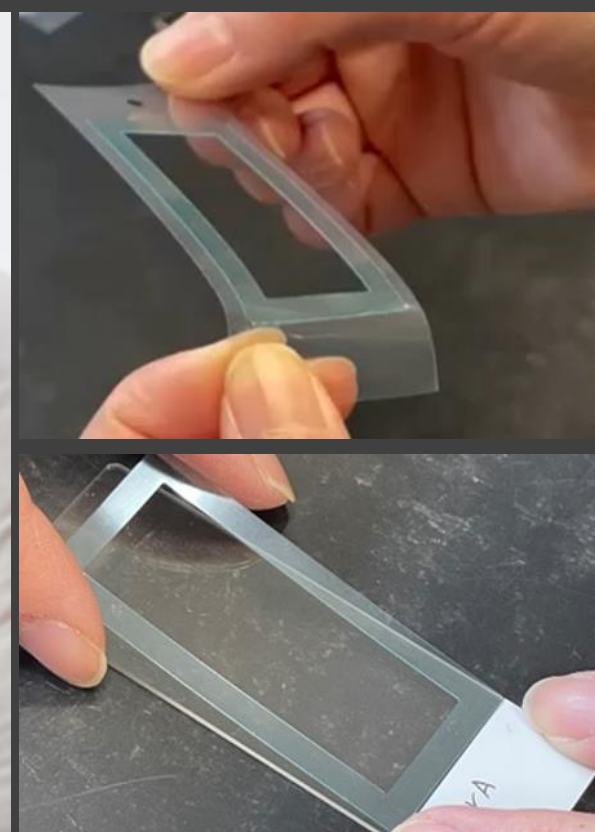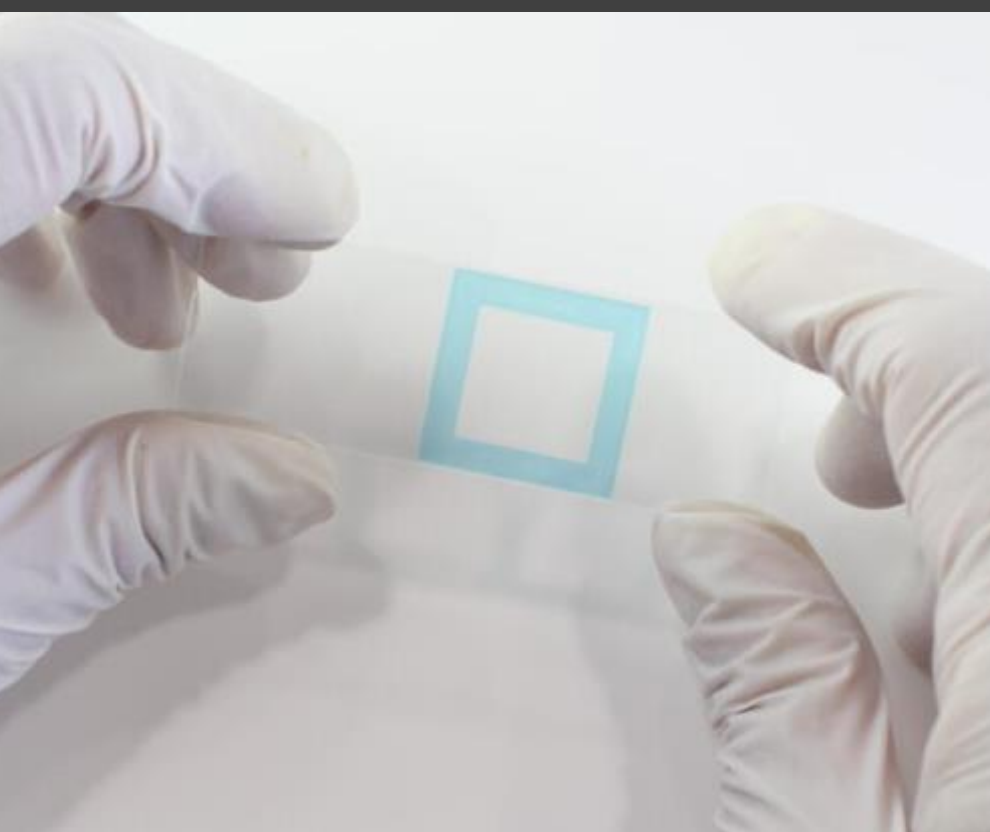

- **Sticker format enclosure for slide-mounted tissue**
- **Water-repellent seal prevents leakage and outflow of antibody or buffer solution in immunostaining**

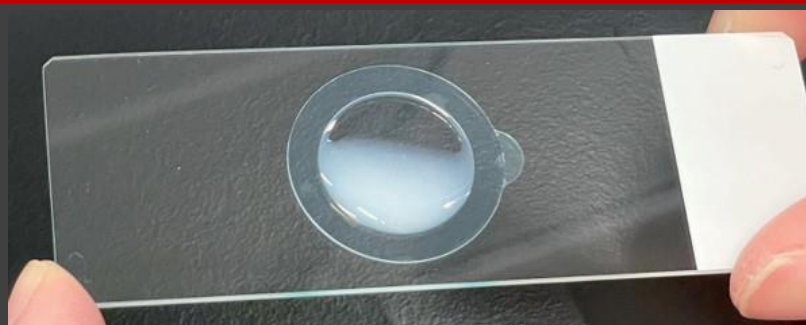


Easy reposition and peel off by peel tab

No need to wipe off moisture before placing the barrier seal

Retains adhesiveness even at high temperatures (70 °C)

6 types are available

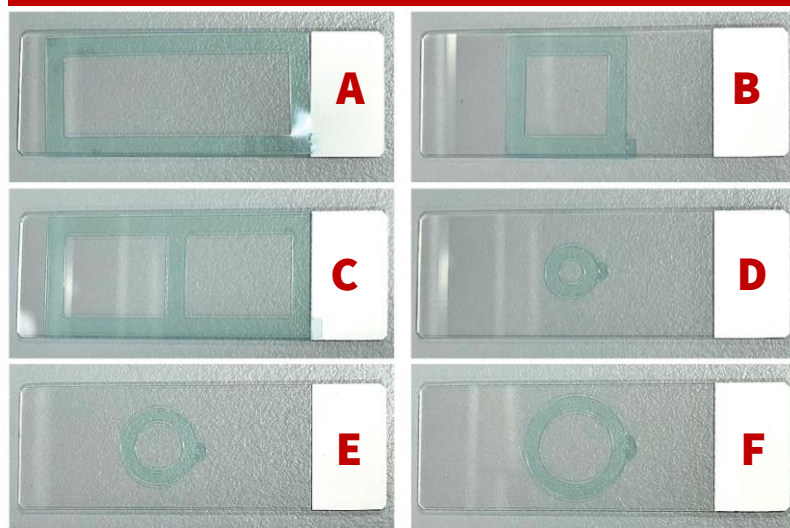

Material: PET (removable adhesive)


***If there are air bubbles under the seal, please push it out.**

How To Use

Section tissue and place on slides



Put on a barrier seal

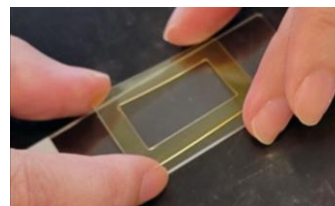

Fixation and blocking




Add primary antibody



Wash and add secondary antibody



Wash, dehydrate and mount

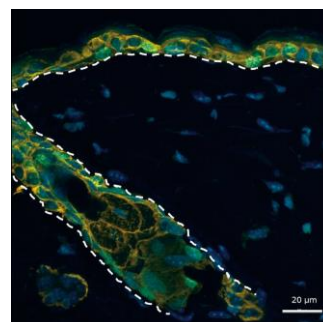

Catalog #	Type	Dimension (mm)		Size
R-CSC091-01	A		Outer: 55×25, Inner: 47×17, Height: 0.155 mm	60 pieces
R-CSC092-01	B		Outer: 25×25, Inner: 17×17, Height: 0.155 mm	90 pieces
R-CSC093-01	C		Outer: 55×25, Inner: 21.5×17, Height: 0.155 mm	60 pieces
R-CSC094-01	D		Outer φ: 11, Inner φ: 5, Height: 0.155 mm	90 pieces
R-CSC095-01	E		Outer φ: 16, Inner φ: 10, Height: 0.155 mm	90 pieces
R-CSC096-01	F		Outer φ: 21, Inner φ: 15, Height: 0.155 mm	90 pieces

Your Local Distributor

Funakoshi Co., Ltd. (JAPAN)

Phone: +81-3-5684-6296

Mail: export@funakoshi.co.jp



\Contact/ \Newsletter Signup/

CSC-2304-F22 (2023.04)